

How to turn off the Cs source

- 1) Turn electrons OFF
- 2) Then turn the electron gun filament to zero (**e- fil.v** button on lower left of keyboard – **red arrow** - and use the **Zoom/Source wheel** – **green arrow**). This may be done quickly.



- 3) To shut the Cs source down:
 - a) go to COLUMNS
 - b) select Stop Source (blue button on right of window)
 - c) select button Cs
 - d) click Apply. Answer OK to the 'log will be updated' message.
- 4) Cs source shut down will proceed automatically and take ~10 minutes. You may leave while this is going on.

